

NOWPak® Bottle Top Adapter Kit

Model No. ND-LD-01 · EMD Part No. 693001-2

Application

This kit is used to connect Optifix® dispensers to NOWPak® containers. Proper installation insures sealed, airtight dispensing.

Instructions For Use

1. Remove air vent cap from Optifix® dispenser. If after reading these instructions you have any questions about assembly or use please contact EMD Chemicals Technical Services for help at 800-222-0342.
2. Install small o-ring on air vent cap and thread back into Optifix® dispenser.
3. Cover air vent hole on cap with sticker to seal NOWPak® liner during dispensing.
4. Install large o-ring onto collar of Optifix® dispenser body.
5. Install PTFE adapter through clearance hole in adapter cap.
6. Remove cap from NOWPak® container.
7. Tighten on adapter cap subassembly by hand.
8. Push flared end of diptube onto Optifix® dispenser valve body nipple.
9. Insert Optifix® dispenser with diptube through connector and lock components together per dispenser instructions.

Note: For Optifix® dispensers 100 ml and larger, tubing must be replaced with larger diameter tubing and the PTFE Adapter modified.

For any application questions call ATMI (800-966-6698).

For questions about the Optifix® dispenser contact EMD Chemicals.

- 1 Optifix® Dispenser (not provided)
- 2 o-ring Seal
- 3 Adapter Cap
- 4 PTFE Adapter
- 5 20L Diptube



EMD Chemicals Inc. ("EMD") warrants that these products are free from defects in material and workmanship for a period of one (1) year (ninety (90) days for seals and gaskets) from the date of delivery. EMD's obligation under this warranty shall be limited to the repair or replacement, at EMD's expense, of products or parts which are proven to EMD's satisfaction to be defective. This warranty does not cover damage due to accident, misuse, negligence, or abnormal use or service. Any alteration or removal of the serial number on these products shall void this warranty. When making warranty claims, the purchaser must, if requested by EMD, provide to EMD the affected product/part with its serial number unaltered and visible. THIS LIMITED WARRANTY SHALL BE IN LIEU OF ALL WARRANTIES OF FITNESS AND MERCHANTABILITY. EMD SHALL NOT BE LIABLE FOR PRODUCTION LOSSES, LOSS OF PROFITS, DIRECT, INDIRECT, INCIDENTAL, SPECIAL, CONSEQUENTIAL OR OTHER DAMAGES OF ANY KIND NO MATTER HOW ARISING.

Prices and specifications are subject to change without notice. We reserve the right to discontinue items without prior notice.

Optifix® is a registered trademark of Walter Graf u. Co, GmbH & Co.

NOWPak® is a registered trademark of ATMI

EMD Chemicals Inc.
480 S. Democrat Road
Gibbstown, NJ 08027
Phone 800-222-0342
Fax 856-423-4389
www.emdchemicals.com
LT 011059 REV 08/05

