



Product Information

Blue Dextran

D5751

Molecular Weight: 2,000,000

CAS Number: 87915-38-6

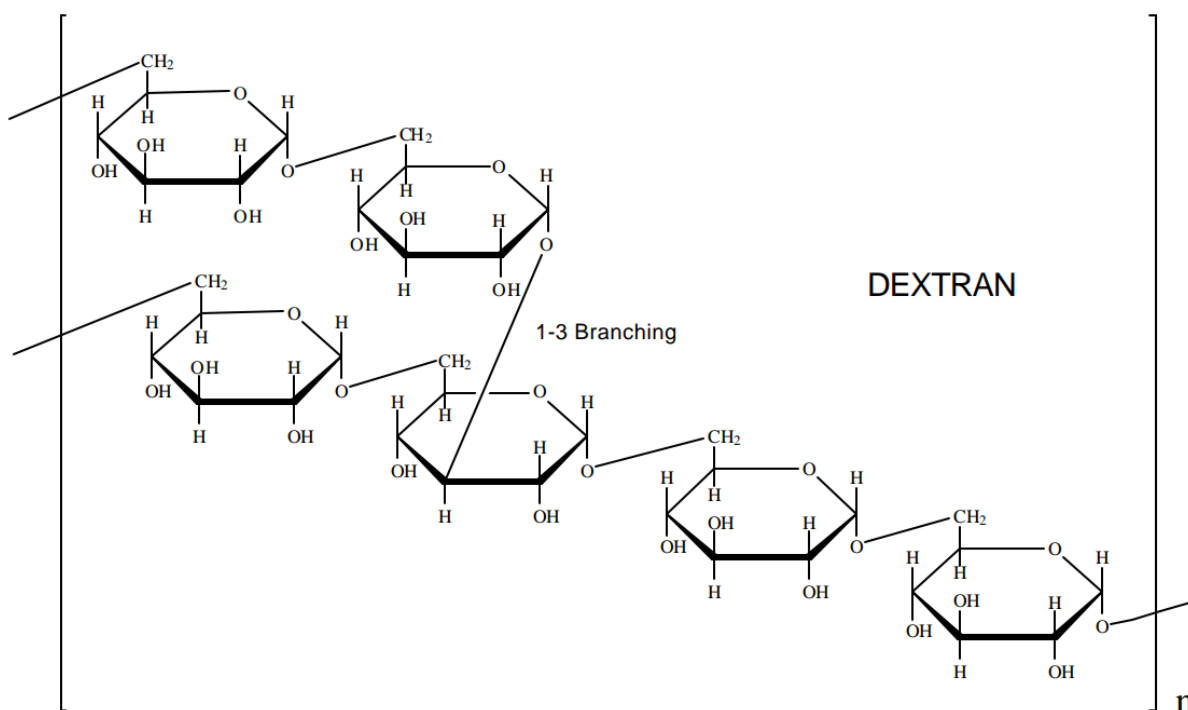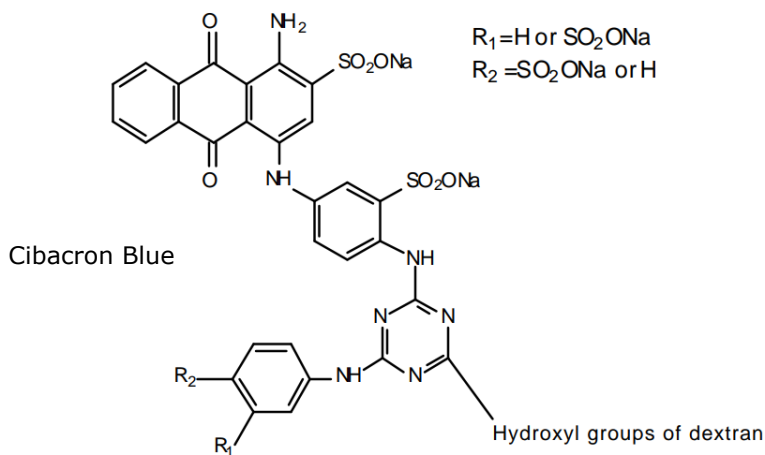

Physical Properties

Absorbance

Lambda max. at approx. 620 nm, E 1% ranged from 8.92 to 9.48

Lambda max. at approx. 380 nm, E 1% ranged from 4.32 to 4.73

Appearance

Dark Blue Powder

Structure

Dextran is a polymer of anhydroglucose. It is composed of approximately 95% alpha-D-(166) linkages. The remaining alpha(163) linkages account for the branching of dextran.^{1,2,3} Conflicting data on the branch lengths implies that the average branch length is less than three glucose units.^{4,5} However, other methods indicate branches of greater than 50 glucose units exist.^{6,7} Native dextran has been found to have a molecular weight (MW) in the range of 9 million to 500 million.^{8,9,10} Dextran with MW greater than 10,000 behave as if they are highly branched. As the MW increases, dextran molecules attain greater symmetry.^{7,11,12} The MW of dextran is measured by one or more of the following methods: low angle laser light scattering (LALLS),¹³ size exclusion chromatography,¹⁴ copper-complexation¹⁵ and anthrone reagent¹⁶ colorimetric reducing-end sugar determination and viscosity.¹¹ The Cibracon Blue Dye of blue dextran is attached randomly to hydroxyl groups along the dextran chain.

Specific Rotation for Dextran

$[\alpha] = +199E^{17}$

Method of Preparation

Our dextrans are derived from *Leuconostoc mesenteroides*, strain B 512. Various MWs are produced by limited hydrolysis and fractionation. Our supplier's exact methods are held proprietary. Fractionation can be accomplished by size exclusion chromatography¹⁴ or ethanol fractionation in which the largest MW dextrans precipitate first.¹⁸ The reactive blue dye is coupled directly to the hydroxyl groups of dextran by a method similar to that used to prepare reactive dye affinity matrices.¹⁹

Stability/Storage as Supplied

This product is given a shelf-life of five years from the date of manufacture.

Solubility/Solution Stability

We test the solubility of Blue Dextran at 50 mg/mL in water. Dextran can be hydrolyzed by strong acids at high temperatures. The terminal reducing end group of dextran can be oxidized in alkaline solutions.¹⁷ Solutions should be prepared fresh whenever possible. Dextran is susceptible to microbial degradation. Solutions should be stored refrigerated.

Applications

The most common use for Blue Dextran is in the determination of the void volume in gel filtration chromatography columns. It is typically used at 1 mg/mL to 2 mg/mL in a buffer such as 50 mM Tris-HCl, pH 7.5 containing 100 mM KCl. The volume of the Blue Dextran sample should be 1% to no more than 3% of the column's bed volume. The shape of the Blue Dextran band as it elutes through the column can also be used to assess the column packing and resolution.

Blue Dextran has also been used as a spectrophotometric substrate for dextranase. Dextranase activity is measured by the increase in absorbance at 610 nm.¹⁹

The affinity of the Cibracon Blue Dye in Blue Dextran for certain proteins has been exploited in protein purification methods. When bound to a protein the Blue Dextran-protein complex will be eluted in the void volume of a Sephadex[®] gel filtration column. This allows the bound protein to be easily separated from other proteins of similar molecular weight. The binding of Blue Dextrans to proteins is dependent on pH and ionic strength.^{20,21}

References

1. Rankin, J.C. and Jeanes, A., J. Am. Chem. Soc., 76, 4435 (1954).
2. Dimler, R.J. et al., J. Am. Chem. Soc., 77, 6568 (1955).
3. Van Cleve, J.W. et al., J. Am. Chem. Soc., 78, 4435 (1956).
4. Lindberg, B. and Svensson, S., Acta. Chem. Scand., 22, 1907 (1968).
5. Larm, O. et al., Carbohydr. Res. 20, 39 (1971).
6. Bovey, F.A., J. Polym. Sci., 35, 167 (1959).
7. Senti, R.F. et al., J. Polym. Sci., 17, 527 (1955).
8. Arond, L.H. and Frank, H.P., J. Phys. Chem., 58, 953 (1954).
9. Elias, H.G., Makromol. Chem., 33, 166 (1959).
10. Antonini, E. et al., Biopolymers, 2, 27 (1964).
11. Granath, K.A., J. Colloid Sci., 13, 308 (1958).
12. Wales, M. et al., J. Polym. Sci., 66, 101 (1979).
13. Allen, P.W., Techniques of Polymer Characterization, Butterworths Scientific Publications, p. 131 (1959).
14. Granath, K.A. and Flodin, P., Makromol. Chem., 48, 160 (1961).
15. Journal of Research of the National Bureau of Standards, 50, 81 (1953).
16. Jermyn., M.A., Anal. Biochem., 68, 332 (1975).

Supplier's Data

17. Ingelman, B. and Halling, M.S., Ark. Kemi., 1, 61 (1949).
18. Kroviarski, Y., J. Chromatog., 449, 403 (1988).
19. Blume, K.G., Biochim. Biophys. Acta, 227, 364 (1971).
20. Swart, A.C.W. and Hemker, H.C., Biochim. Biophys. Acta, 222, 692 (1970).

Notice

We provide information and advice to our customers on application technologies and regulatory matters to the best of our knowledge and ability, but without obligation or liability. Existing laws and regulations are to be observed in all cases by our customers. This also applies in respect to any rights of third parties. Our information and advice do not relieve our customers of their own responsibility for checking the suitability of our products for the envisaged purpose.

The information in this document is subject to change without notice and should not be construed as a commitment by the manufacturing or selling entity, or an affiliate. We assume no responsibility for any errors that may appear in this document.

Technical Assistance

Visit the tech service page at SigmaAldrich.com/techservice.

Terms and Conditions of Sale

Warranty, use restrictions, and other conditions of sale may be found at SigmaAldrich.com/terms.

Contact Information

For the location of the office nearest you, go to SigmaAldrich.com/offices.

The life science business of Merck KGaA, Darmstadt, Germany operates as MilliporeSigma in the U.S. and Canada.

MilliporeSigma, and Sigma-Aldrich are trademarks of Merck KGaA, Darmstadt, Germany or its affiliates. All other trademarks are the property of their respective owners. Detailed information on trademarks is available via publicly accessible resources.

© 1997-2024 Merck KGaA, Darmstadt, Germany and/or its affiliates. All Rights Reserved.
D5751pis Rev 11/24

**MILLIPORE
SIGMA**