

Product Information

β-Glucuronidase from *Escherichia coli*, recombinant from overexpressing *Escherichia coli* BL21

G8295

Storage Temperature -20 °C

CAS RN 9001-45-0

EC 3.2.1.31

Synonyms: β-D-Glucuronide glucuronosohydrolase; GUS

Product Description

Glucuronidation, conjugation with glucuronic acid, by the human UDP-glucuronosyltransferase (UGT) family of enzymes plays an important role in the metabolic fate of many drugs and other xenobiotics. This biosynthetic reaction also has a role in the conjugation and excretion of endogenous substrates, such as steroids, bilirubin, and bile acids.¹ UGT activity results in the conjugation of glucuronic acid to substrates containing sulfhydryl, hydroxyl, aromatic amino, or carboxylic acid moieties. The glucuronides formed are more polar (water soluble) than the parent organic substrate and are generally excreted through the kidney.

β-glucuronidase catalyzes the reaction:

$$\beta\text{-D-glucuronoside} + \text{H}_2\text{O} \leftrightarrow \text{an alcohol} + \text{D-glucuronate}$$

β-Glucuronidase from *E. coli* is used for the enzymatic hydrolysis of β-glucuronides in urine and other fluids. It does not hydrolyze α-glucuronides or β-glucosides.² The enzyme from *E. coli* has a high rate of hydrolytic activity and it retains this activity during hydrolysis better than similar enzymes that are more sensitive to changes in the concentration of β-glucuronide conjugates. The enzyme preparation from *E. coli* has been shown to be useful for determining the presence of androsterone, 17-hydroxycorticosteroids, and estriol in urine.³ The optimal conditions for the enzymatic hydrolysis of α-hydroxytriazolam, one of the major metabolites of triazolam in human urine, were determined using β-glucuronidase from *E. coli*. It was found that a 90 minute incubation of 1 mL of urine with 100 units of the enzyme at 37 °C and pH 5.5-7.8, effectively

hydrolyzed the α-hydroxytriazolam given at the clinical dose.⁴

This β-Glucuronidase product from *E. coli* is supplied as a powder lyophilized from 10 mM potassium phosphate. Sucrose is added as a stabilizer.

Molecular Weight: ~290 kDa (tetramer)⁵
68,259 Da (monomer)⁶

Optimal pH⁵: 6-7

Isoelectric point (pI)⁵: 4.8

Inhibitors: D-glucuronic acid (Cat. No. G5269)
D-galacturonic acid
D-glucaro-1,4-lactone

Substrates:

5-Bromo-6-chloro-3-indolyl β-D-glucuronide	B4532
6-Bromo-2-naphthyl β-D-glucuronide	B6519
5-Bromo-4-chloro-3-indolyl β-D-glucuronide sodium salt tablet	B8174
8-Hydroxyquinoline glucuronide	H1254
4-Methylumbelliferyl β-D-glucuronide	M5664
4-Nitrophenyl β-D-glucuronide	73677

Glucuronidase Activity: ≥ 10,000,000 units/gm protein

Unit Definition: One Sigma or modified Fishman unit will liberate 1.0 µg of phenolphthalein from phenolphthalein glucuronide per hour at 37 °C at pH 6.8 (30 minute assay).

Unlike the enzyme preparation from snail (*Helix pomatia*) that naturally contains β-glucuronidase and sulfatase activities in almost equal amounts, the preparation of β-glucuronidase from *E. coli* is essentially free of sulfatase activity.

Precautions and Disclaimer

For R&D use only. Not for drug, household, or other uses. Please consult the Safety Data Sheet for information regarding hazards and safe handling practices.

Solubility

When reconstituted to 5 mg/ml in 75 mM phosphate buffer, pH 6.8, a clear to slightly hazy solution results. Regardless of the cloudiness, the enzyme is active and should be useable for metabolite hydrolysis.

Storage/Stability

The product as supplied should be stored at -20 °C. A solution in 75 mM phosphate buffer, pH 6.8, (≥5 mg/ml) may be stored at -20 °C for up to 2 months with little or no loss of activity.

References

1. Tephly, T.R., et al., *Adv. Pharmacol.*, **42**, 343-346 (1998).
2. *Methods of Enzymatic Analysis*, vol. 2, Hans Ulrich Bergmeyer, Ed., Academic Press, NY, 460-461, 929-943 (1974).
3. Graef, V., et al., *Clin. Chem.*, **23**, 532-535 (1977).
4. Tsujikawa, K., et al., *J. Health Sci.*, **50**, 286-289 (2004).
5. Kim, D-H, et al., *Biol. Pharm. Bull.*, **18**, 1184-1188 (1995).
6. Jefferson, R.A., et al., *Proc. Natl. Acad. Sci.*, **83**, 8447-8451 (1986).

AC, RBG, MAM, KKK 01/09-1

Notice

We provide information and advice to our customers on application technologies and regulatory matters to the best of our knowledge and ability, but without obligation or liability. Existing laws and regulations are to be observed in all cases by our customers. This also applies in respect to any rights of third parties. Our information and advice do not relieve our customers of their own responsibility for checking the suitability of our products for the envisaged purpose.

The information in this document is subject to change without notice and should not be construed as a commitment by the manufacturing or selling entity, or an affiliate. We assume no responsibility for any errors that may appear in this document.

Technical Assistance

Visit the tech service page at SigmaAldrich.com/techservice.

Terms and Conditions of Sale

Warranty, use restrictions, and other conditions of sale may be found at SigmaAldrich.com/terms.

Contact Information

For the location of the office nearest you, go to SigmaAldrich.com/offices.

The life science business of Merck KGaA, Darmstadt, Germany operates as MilliporeSigma in the U.S. and Canada.

MilliporeSigma, and Sigma-Aldrich are trademarks of Merck KGaA, Darmstadt, Germany or its affiliates. All other trademarks are the property of their respective owners. Detailed information on trademarks is available via publicly accessible resources.

© 2023 Merck KGaA, Darmstadt, Germany and/or its affiliates. All Rights Reserved.

G8295PIS Rev 02/23