usage in vitro

Anaerocult® T

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
16820 Microbiology

Anaerocult® T

Contents 25 Anaerocult® T
25 Special incubation bags

Application Production of an anaerobic environment in agar slant tubes for the cultivation and transport of Obligate and facultative anaerobes.


Principle Anaerocult® T contains components that chemically bind any oxygen present within a short space of time and also release carbon dioxide thus creating an anaerobic atmosphere.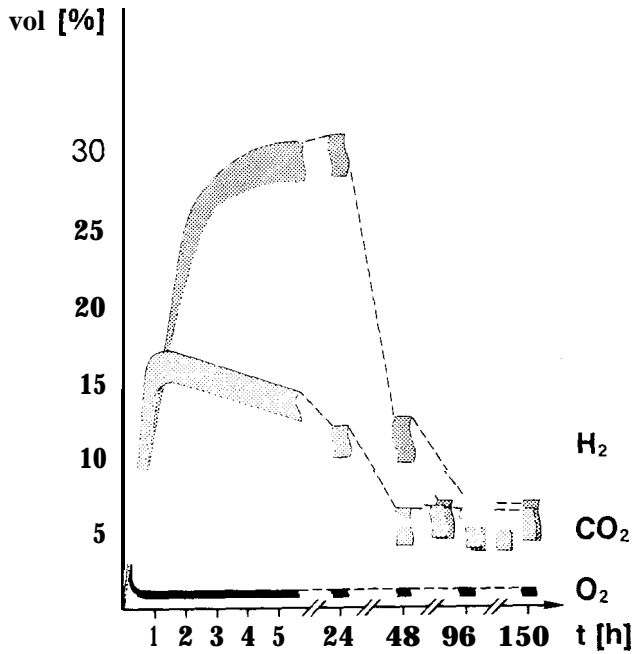


Composition Kieselguhr
Iron powder
Citric acid
Sodium carbonate

Directions Anaerocult® T is placed in the special incubation bag together with the agar slant tube and the anaerobiosis indicator Anaerostest® (Cat. No. 15112).


Procedure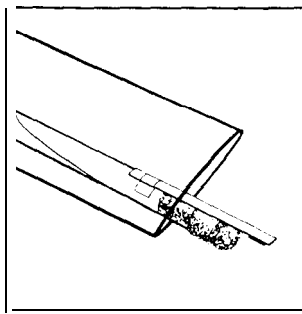


1 Moisten the reaction zone of Anaerostest™ with water and stick the anaerobiosis indicator on to one slant tube (the reaction zone must hang freely in the open space) (Fig. 1). Place 1-3 agar slant tubes (sealed with cotton wool plugs) in the special incubation bag (Fig. 1).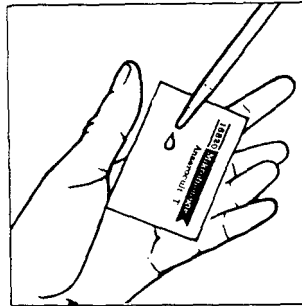

2 Moisten Anaerocult® T with 3 ml water (Fig. 2).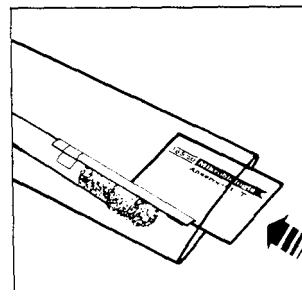

3 Place the moistened Anaerocult® T immediately in the special incubation bag (Fig. 3).

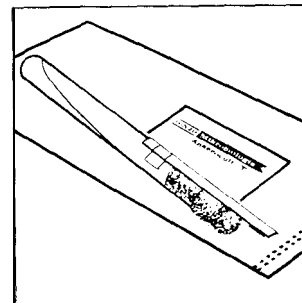4 Seal the special incubation bag with a foil sealing device (it is advisable to make 2 seals)* (Fig. 4).

Stability See expiry date.

Storage Seal tightly and protect from moisture (seal the plastic bag well after removing Anaerocult® T). Recommended storage temperature: +15 °C ↔ +25 °C

* Anaerobiosis is indicated by the colour change of the Anaerostest®-strip from blue to white after about 4 hours.

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