

Product Information

Protease Inhibitor Cocktail

For general use

P2714

Product Description

Crude cell extracts contain various endogenous enzymes, such as proteases and phosphatases, which can degrade proteins in the extracts. The best way to increase the yield of intact proteins is to add inhibitors of those enzymes known to be present.

The P2714 protease inhibitor cocktail has been optimized and tested for general use. P2714 is a mixture of water-soluble protease inhibitors with a broad specificity for the inhibition of serine, cysteine, and metalloproteases.

P2714 is supplied as a lyophilized powder. The inhibitors in P2714 are as follows, with respective specific inhibitor targets and target classes of each inhibitor listed:

- AEBSF [4-(2-Aminoethyl)benzenesulfonyl fluoride hydrochloride]: serine proteases, such as trypsin, chymotrypsin, plasmin, kallikrein and thrombin
- Aprotinin: serine proteases, such as trypsin, chymotrypsin, plasmin, and kallikrein; human leukocyte elastase, but not pancreatic elastase
- Bestatin hydrochloride: aminopeptidases, such as leucine aminopeptidase and alanyl aminopeptidase¹⁻⁴
- E-64 [*N*-(trans-Epoxy succinyl)-L-leucine 4-guanidinobutylamide]: cysteine proteases, such as calpain, papain, cathepsin B, and cathepsin L
- EDTA: metalloproteases
- Leupeptin hemisulfate salt: serine proteases and cysteine proteases, such as plasmin, trypsin, papain, and cathepsin B

Several theses⁵ and dissertations⁶⁻²⁰ have cited use of product P2714 in their protocols.

Storage/Stability

Store the lyophilized powder at -20 °C. The lyophilized powder is stable for 4 years when stored at -20 °C.

Precautions and Disclaimer

For R&D use only. Not for drug, household, or other uses. Please consult the Safety Data Sheet for information regarding hazards and safe handling practices.

Preparation Instructions

One bottle prepares 100 mL of cocktail solution.

The product is packaged in a 10 mL serum vial with a crimp cap. Dissolve the contents of the vial in 10 mL of working buffer, then transfer to another container for dilution to 100 mL.

Recommended Usage

- One mL of the P2714 cocktail solution is recommended for the inhibition of endogenous enzymes equivalent to 1 mg of USP pancreatin.
- One bottle is recommended for the inhibition of proteases present in a maximum of 20 g of cell extract.
- **Note:** Not all lysates contain the same levels of endogenous enzymes. It may be necessary to adjust the volume of cocktail required.

References

1. Umezawa, H., *Ann. Rev. Microbiol.*, **36**, 75-99 (1982).
2. Aoyagi, T. *et al.*, *Biochem. Int.*, **9(4)**, 405-411 (1984).
3. Aoyagi, T., and Umezawa, H., *Acta Biol. Med. Ger.*, **40(10-11)**, 1523-1529 (1981).
4. Mumford, R.A. *et al.*, *Biochem. Biophys. Res. Comm.*, **103(2)**, 565-572 (1981).
5. Rennie, Keith, "The Role of DAP-Kinase in Modulating Vascular Endothelial Cell Function Under Fluid Shear Stress". Purdue University, M.S. thesis, p. 26 (2010).

6. Mimeault, Caroline, "Human Pharmaceuticals in the Environment: Effects of Fibrate Drugs on Two Fish Species". University of Ottawa, Ph.D. dissertation, p. 44 (2008).
7. Shulla, Ana, "Cofactors in Coronavirus Entry". Loyola University Chicago, Ph.D. dissertation, pp. 49, 50 (2011).
8. Yoon, Jennifer Ruth, "The Role of Kinesin Motor Proteins in Mammary Epithelial And Breast Tumor Cell Microtentacles". University of Maryland Baltimore, Ph.D. dissertation, pp. 20, 89 (2011).
9. Gibrel, Gehan G.F., "The Role of ADAMTS-1, -4 and -5 in Multiple Sclerosis". Sheffield Hallam University, Ph.D. dissertation, p. 86 (2012).
10. Winsor, Tiffany Sabin, "A Novel DNA Damage Response Network Associated with the CTD of RNA Polymerase II". Duke University, Ph.D. dissertation, p. 64 (2012).
11. Broadwater, Laurie A., "Differential Proteomic Investigations of Normal Appearing Gray Matter in Multiple Sclerosis and Control Post-Mortem Brain Tissue". Kent State University, Ph.D. dissertation, p. 36 (2013).
12. Bjerke, Maureen Ann, "Mechanosensitive adhesive networks guide collective migration of *Xenopus* mesendoderm". University of Virginia, Ph.D. dissertation, p. 129 (2014).
13. McGuiness, Graham Ripley, "Interleukin-6 and Exercise-Induced Cardioprotection". Auburn University, Ph.D. dissertation, pp. 29, 77 (2014).
14. Bedoya, Flor Edith Acevedo, "The wear and tear of feeding: induced plant defenses by fall armyworm herbivory and counter insect adaptation mechanisms". The Pennsylvania State University, Ph.D. dissertation, pp. 50, 51, 127 (2016).
15. Ko, Dae Kwan, "Clock-regulatory networks contributes to growth vigor in maize hybrids". University of Texas at Austin, Ph.D. dissertation, pp. 93, 96 (2016).
16. Figueirinhas do Amaral, Maria Clara, "Geographical Variation of Freeze Tolerance in the Wood Frog, *Rana sylvatica*: The Role of Hepatic Glycogen Metabolism". Miami University, Ph.D. dissertation, pp. 12, 42 (2017).
17. Santiago-Frangos, Andrew, "Modular disordered domains autoregulate the RNA binding and annealing activities of the bacterial RNA chaperone Hfq". Johns Hopkins University, Ph.D. dissertation, pp. 159, 181, 182 (2018).
18. Piszkiwicz, Samantha, "Tardigrade Disordered Proteins as Potential Excipients for Biologics". University of North Carolina at Chapel Hill, Ph.D. dissertation, pp. 34, 63 (2019).
19. Schlütermann, David, "Autophagy: Molecular insights into its role and therapeutic potential in bladder cancer and neurodegeneration". Heinrich-Heine-Universität Düsseldorf, Dr. rer. nat. dissertation, pp. 44, 111 (2020).
20. Kunselman, Jennifer M., "Regulation of Opioid Receptor Trafficking and Signaling by Opioid Peptides ". University of Michigan, Ph.D. dissertation, pp. 43, 44, 46, 47, 51 (2021).

Notice

We provide information and advice to our customers on application technologies and regulatory matters to the best of our knowledge and ability, but without obligation or liability. Existing laws and regulations are to be observed in all cases by our customers. This also applies in respect to any rights of third parties. Our information and advice do not relieve our customers of their own responsibility for checking the suitability of our products for the envisaged purpose.

The information in this document is subject to change without notice and should not be construed as a commitment by the manufacturing or selling entity, or an affiliate. We assume no responsibility for any errors that may appear in this document.

Technical Assistance

Visit the tech service page at [SigmaAldrich.com/techservice](https://www.sigmaaldrich.com/techservice).

Standard Warranty

The applicable warranty for the products listed in this publication may be found at [SigmaAldrich.com/terms](https://www.sigmaaldrich.com/terms).

Contact Information

For the location of the office nearest you, go to [SigmaAldrich.com/offices](https://www.sigmaaldrich.com/offices).

The life science business of Merck KGaA, Darmstadt, Germany operates as MilliporeSigma in the U.S. and Canada.

MilliporeSigma, and Sigma-Aldrich are trademarks of Merck KGaA, Darmstadt, Germany or its affiliates. All other trademarks are the property of their respective owners. Detailed information on trademarks is available via publicly accessible resources.

© 2022 Merck KGaA, Darmstadt, Germany and/or its affiliates. All Rights Reserved.

P2714dat Rev 06/22 EM,PHC,GCY

**MILLIPORE
SIGMA**