



3050 Spruce Street
Saint Louis, Missouri 63103 USA
Telephone (800) 325-5832 (314) 771-5765
Fax (314) 286-7828
email: techserv@sial.com
sigma-aldrich.com

Product Information

Ubiquitin Activating Enzyme (E1) from rabbit

Product Number **U1758**

Storage Temperature $-70\text{ }^{\circ}\text{C}$

Product Description

Ubiquitin activating enzyme (E1) is a 110 kDa protein isolated from rabbit.

The ubiquitin proteolytic system plays an important role in a broad array of basic cellular processes. Among these are regulation of cell cycle, modulation of the immune and inflammatory responses, development and differentiation, and control of signal transduction pathways. These complex processes are controlled via specific degradation of a single or a subset of proteins. Degradation of a protein by the ubiquitin system involves two successive steps, conjugation with multiple moieties of ubiquitin and degradation of the ubiquitin tagged protein by the 26S proteasome.¹

Ubiquitin activating enzyme (E1) performs the initial step in the formation of the ubiquitin-protein isopeptide bond. E1 catalyzes the activation of the C-terminal carboxyl group of ubiquitin by forming a high energy thioester bond in an ATP-dependent manner.² This activated ubiquitin is then transferred to a lysine of the target protein via the E2/E3 conjugation cascade. It is the critical step for the initiation of any *in vitro* conjugation reactions.³ A working concentration of 50 to 200 nM is recommended to support *in vitro* conjugation.

This product is supplied as a solution in 50 mM HEPES, pH 7.6.

Purity: minimum 95% (SDS-PAGE)

Precautions and Disclaimer

This product is for laboratory research use only. Please consult the Material Safety Data Sheet for information regarding hazards and safe handling practices.

Storage/Stability

This product ships on dry ice and storage at $-70\text{ }^{\circ}\text{C}$ is recommended. The product is stable through multiple freeze/thaw cycles.

References

1. Ciechanover, A., The Ubiquitin-mediated proteolytic pathway: mechanisms of action and cellular physiology. *Biol. Chem. Hoppe-Seyler*, **375**, 565-581 (1994).
2. Ciechanover, A., et al., "Covalent affinity" purification of ubiquitin-activating enzyme. *J. Biol. Chem.*, **257**, 2537-2542 (1982).
3. Haas, A., et al., Ubiquitin-activating enzyme. Mechanism and role in protein-ubiquitin conjugation. *J. Biol. Chem.*, **257**, 2543-2548 (1982).

Manufactured for Sigma by Boston Biochem., Inc.

PMH/MAM 9/02

Sigma brand products are sold through Sigma-Aldrich, Inc.

Sigma-Aldrich, Inc. warrants that its products conform to the information contained in this and other Sigma-Aldrich publications. Purchaser must determine the suitability of the product(s) for their particular use. Additional terms and conditions may apply. Please see reverse side of the invoice or packing slip.