

93745 TSC Agar (Tryptose Sulfite Cycloserine Agar, Perfringens Agar)

For the isolation and enumeration of vegetative forms as well as spores from *Clostridium perfringens* in food, clinical specimens and other material acc. to Harmon, et al. (1971).

Composition:

Ingredients	Grams/Litre
Meat peptone	15.0
Soya peptone	5.0
Yeast extract	5.0
Ammonium ferric citrate	1.0
Sodium disulfite	1.0
Agar	15.0
Final pH 7.6 +/- 0.2 at 25°C	

Store prepared media below 8°C, protected from direct light. Store dehydrated powder, in a dry place, in tightly-sealed containers at 2-25°C.

Appearance: Faintly beige, homogeneous, free flowing powder.
Gelling: Firm.
Colour and Clarity: Light brownish-yellow coloured, clear gel forms in petri plates.

Directions:

Base Agar:

Dissolve 42 g in 1 litre distilled water. Sterilize by autoclaving at 121°C for 15 minutes. Cooling to approx. 50°C.

TSC Agar:

To 1 litre Base Agar add 400 mg rehydrated D-cycloserine (Sigma 30020) as a sterile filtered aqueous solution or 2 vials of Perfringens T.S.C. Supplement (P9352) and 50 ml Egg Yolk Emulsion (17148). Mix well and pour into petri plates.

TSC Agar egg yolk free (also used as overlay agar):

Add 400 mg rehydrated D-cycloserine (Sigma 30020) or 2 vials of Perfringens T.S.C. Supplement (P9352) to 1 litre Base Agar and mix well.

TSC Fluorescence Agar:

Add to 1 litre Base Agar 5.5 g rehydrated TSC Agar supplement (80548) as a sterile filtered aqueous solution. Mix well and pour plates.

SFP Agar (Shahidi-Ferguson Perfringens Agar):

To 1 litre Base Agar 2 vials of Perfringens S.F.P. Supplement (P9477) and 50 ml Egg Yolk Emulsion (17148) are added. Mix well and pour into sterile petri plates.



Principle and Interpretation:

Meat peptone, soya peptone and yeast extract provide essential nutrients and vitamins for the development of clostridia. The H₂S-positive bacteria reduce the sulfite (from Sodium disulfite) in the culture medium to sulfide, which forms a black salt with ammonium ferric citrate (FeS).

Egg yolk promotes the growth of *Cl. perfringens* but for counting of plates it can be an advantage to add no egg yolk and have smaller colonies [4,5,6].

No lecithinase activity can be detected without egg yolk but not all *Clostridium perfringens* strains show an opaque zone after overnight incubation [4]. So, both black lecithinase-positive and black lecithinase-negative colonies should be considered as presumptive *Clostridium perfringens* on TSC or SFP Agar and need further confirmatory tests like nitrate reduction, lactose fermentation, gelatin liquefaction and the absence of motility [8].

Cycloserine inhibits the accompanying bacterial flora. It is the reason why some inhibited colonies appear smaller. It also reduces the diffuse and disturbing blackening around the *Cl. perfringens* colonies. All this can simplify the counting of plates with high numbers of colonies. Higher counts have been demonstrated by using a pour plate technique. The differences are thought to be due to exposure of the *Cl. perfringens* cells to high oxygen tension in the surface plating procedure.

TSC Agar Supplement contains Sucrose, D-cycloserine and 4-Methylumbelliferylphosphate (MUP). *Clostridium perfringens* possess the enzyme phosphatase, which cleaves MUP and results in setting free the fluorogen 4-methylumbelliferone. The fluorogen can be detected under a long wavelength UV lamp (360 nm). Sucrose is a fermentable carbohydrate source for many clostridia. It supports the growth and germination of clostridia. With the addition of indicator e.g. 0.04 g/l bromocresol purple it is possible to differentiate further (Sucrose positives: *Cl. perfringens*, *Cl. baratii*, *Cl. paraputrificum*, *Cl. tertium*; Sucrose negatives: *Cl. biferentans*, *Cl. difficile*, *Cl. sporogenes*).

The addition of kanamycin sulfate (12 mg/l; 60615) and polymyxin B sulphate (30,000 IU/l ≈ 3.6 mg/l; 81334) present in the Perfringens S.F.P. Supplement gives a better recovery rate than with cycloserine but still a high degree of selectivity and specificity for *Cl. perfringens* [1,2,3]. It was observed that more unspecific colonies were found on SFP Agar

For incubation are diverse temperatures (35-45°C) recommended. A higher temperature increase the background growth but result in a lower recovery rate[13].

Cultural characteristics after 18-24 hours at 44°C (under anaerobic conditions).

Organisms (ATCC)	Growth	Black colony	Lecithinase activity
<i>Clostridium perfringens</i> (13124)	+++	+	+
<i>Clostridium tetani</i> (19406)	-/++	-	
<i>Clostridium novyi</i> (10543)	-/++	-	
<i>Pseudomonas aeruginosa</i> (27853)	-/+	-	
<i>Bacillus cereus</i> (11778)	-/+	-	



References:

1. S.A. Shahidi, A.R. Ferguson, New quantitative, qualitative and confirmatory media for rapid analysis food for *Clostridium perfringens*, Appl. Microbiol., 21, 500-506 (1971)
2. S.M. Harmon, et al., Appl. Microbiol. 21, 922-927 (1971)
3. S.M. Harmon, et al., Appl. Microbiol. 22, 688-692 (1971)
4. A.H.W. Hauschild, R. Hilsheimer, Evaluation and modifications of media for enumeration of *Clostridium perfringens*, Appl. Microbiol., 27, 78-82 (1974)
5. A.H.W. Hauschild, R. Hilsheimer, Enumeration of food-borne *Clostridium perfringens* in egg yolk-free tryptose-sulfite-cycloserine agar. Appl. Microbiol. 27, 521-526 (1974)
6. A.H.W. Hauschild, R. Hilsheimer, Enumeration of faecal *Clostridium perfringens* in egg yolkfree tryptose-sulphite cycloserine agar. Appl. Microbiol. 27, 527-530 (1974)
7. D.S. Orth, Comparison of sulfite-polymyxin-sulfadiazine medium and tryptose-sulfite-cycloserine medium without egg yolk for recovering *Clostridium perfringens*, Appl. Envir. Microbiol., 33, 986 (1977)
8. A.H.W. Hauschild, R.J. Gilbert, S.M. Harmon, M.F. O'Keefe, R. Vahlfeld, ICMSF methods studies. VIII. Comparative study for the enumeration of *Clostridium perfringens* in foods, Can. J. Microbiol., 23, 884-892 (1977)
9. International Organization for Standardization (ISO), Meat and meat products: Enumeration of *Clostridium perfringens*, Draft ISO/TC 34/SC 6 (1978)
10. T.J. Easterbrook, P.A. West, Isolation and enumeration of sulphite-reducing *Clostridium* spores and Group D faecal streptococci from oysters, J. Appl. Bact. 62, 413 (1987)
11. D.P. Sartory, Membrane filtration enumeration of faecal clostridia and *C. perfringens* in water, Wat. Res. 20, 1255 (1986)
12. American Public Health Association, Compendium of methods for the microbiological examination of foods, 3rd (1992)
13. D.P. Sartory, M. Field, S.M. Curbishley, A.M. Pritchard, Evaluation of two media for the membrane filtration enumeration of *Clostridium perfringens* from water, Appl. Microbiol. 27 (6), 323-327 (1998)

Precautions and Disclaimer

This product is for R&D use only, not for drug, household, or other uses. Please consult the Material Safety Data Sheet for information regarding hazards and safe handling practices.

The vibrant M, Millipore, and Sigma-Aldrich are trademarks of Merck KGaA, Darmstadt, Germany or its affiliates. Detailed information on trademarks is available via publicly accessible resources.
© 2018 Merck KGaA, Darmstadt, Germany and/or its affiliates. All Rights Reserved.

The life science business of Merck KGaA, Darmstadt, Germany operates as MilliporeSigma in the US and Canada.

