



3050 Spruce Street  
 Saint Louis, Missouri 63103 USA  
 Telephone 800-325-5832 • (314) 771-5765  
 Fax (314) 286-7828  
 email: techserv@sial.com  
 sigma-aldrich.com

## Product Information

### Defined Medium Supplements

Product Number	Product Name	Pkg. Size	Reconstitution	Working Range	Storage	References
A 8661	d-ALDOSTERONE	5 mg	To prepare 100 µg/ml stock solution; add 5.0 ml abs. ethanol, gently swirl to dissolve, add 45 ml sterile medium.	2-10 µg/ml	powder-RT stock-frozen in working aliquots, avoid repeated freeze/thaw	1, 2, 10, 54, 55, 56
D 8893	DEXAMETHASONE	1 mg	To prepare 20 µg/ml stock solution; add 1.0 ml abs. ethanol, gently swirl to dissolve, add 49 ml sterile medium.	4-500 ng/ml	powder 2-8°C; stock-frozen in working aliquots, avoid repeated freeze/thaw	11, 47
D 7149	DIHYDRO-TESTOSTERONE	50 mg	To prepare 1 mg/ml stock solution; add 1.0 ml abs. ethanol, gently swirl to dissolve, add 49 ml sterile medium.	2-100 ng/ml	powder-RT stock-frozen in working aliquots, avoid repeated freeze/thaw	
E 2257	β-ESTRADIOL	1 mg	To prepare 20 µg/ml stock solution; add 1.0 ml abs. ethanol, gently swirl to dissolve, add 49 ml sterile medium.	0.2-10 ng/ml	powder-RT stock-frozen in working aliquots, avoid repeated freeze/thaw	12, 34, 36, 42
G 3157	GLUCAGON	2 mg	To prepare 100 µg/ml stock solution; add 20 ml sterile acidified H <sub>2</sub> O (pH≤3.0), swirl to dissolve.	0.03-10 µg/ml	powder -0°C; stock-frozen in working aliquots, avoid repeated freeze/thaw	13
H 0135	HYDROCORTISONE	1 mg	To prepare 50 µg/ml stock solution; add 1.0 ml abs. ethanol, gently swirl to dissolve, add 19 ml sterile medium	0.004-5 µg/ml	powder-RT; stock-frozen in working aliquots, avoid repeated freeze/thaw	8, 14, 35, 36
I 1882	INSULIN	100 mg	To prepare 10 mg/ml stock solution; add 10 ml of acidified H <sub>2</sub> O (pH≤2.0) - prepared by addition of glacial acetic acid (approx. 0.1 ml).	0.001-20 µg/ml	powder -0°C; solution stable 2-8°C, 1 year	3, 4, 15
I 1884	INSULIN-TRANSFERRIN	25 mg, 25 mg,	To prepare a stock solution, dissolve in 5 ml sterile acidified H <sub>2</sub> O, swirl to dissolve,	5 µg, 5 µg,	powder -0°C; stock-frozen in working aliquots, avoid	3, 4

SODIUM SELENITE

25  $\mu$ g

follow with 45 ml sterile water. Stock solution makes 5 L medium final volume.

5 ng/ml

repeated freeze/thaw; solution stable 2-8°C, 30 days

**Defined Medium Supplements** continued

L 8384	LINOLEIC ACID-BOVINE SERUM ALBUMIN (LA/BSA)	500 mg	To prepare 10 mg/ml LA/BSA stock solution, add 50 ml sterile culture medium, gently swirl to dissolve.	1 mg/ml	powder 2-8°C; stock-frozen in working aliquots, avoid repeated freeze/thaw; solution stable 2-8°C, 30 days	3, 27, 30, 53
L 9510	LUTEINIZING HORMONE RELEASING HORMONE	0.1 mg	To prepare 10 µg/ml stock solution; add 10 ml sterile culture medium, gently swirl to dissolve.	0.5-40 ng/ml	powder -0°C; stock-frozen in working aliquots, avoid repeated freeze/thaw	3, 9, 16, 39, 48
P 6149	PROGESTERONE	1.0 mg	To prepare 20 µg/ml stock solution; add 1.0 ml abs. ethanol, gently swirl to dissolve, add 49 ml sterile medium.	0.1-20 ng/ml	powder-RT stock-frozen in working aliquots, avoid repeated freeze/thaw	17, 49
P 6657	PROSTAGLANDIN D2	1.0 mg	To prepare 50 µg/ml stock solution; add 1.0 ml abs. ethanol, gently swirl to dissolve, add 19 ml sterile medium.	0.25-100 ng/ml	powder -0°C; stock-frozen in working aliquots, avoid repeated freeze/thaw	18, 44, 46
P 7527	PROSTAGLANDIN E1	1.0 mg	To prepare 50 µg/ml stock solution; add 1.0 ml abs. ethanol, gently swirl to dissolve, add 19 ml sterile medium.	0.25-100 ng/ml	powder -0°C; stock-frozen in working aliquots, avoid repeated freeze/thaw;	19, 31, 32, 43, 44, 46
P 6532	PROSTAGLANDIN E2	1.0 mg	To prepare 50 µg/ml stock solution; add 1.0 ml abs. ethanol, gently rotate to dissolve, add 19 ml sterile medium	0.25-100 ng/ml	powder -0°C; stock-frozen in working aliquots, avoid repeated freeze/thaw;	18, 44, 46
P 7652	PROSTAGLANDIN F2	1.0 mg	To prepare 50 µg/ml stock solution; add 1.0 ml abs. ethanol, gently swirl to dissolve, add 19 ml sterile medium.	0.25-100 ng/ml	powder -0°C stock-frozen in working aliquots, avoid repeated freeze/thaw; solution stable 2-8°C, 30 days	19, 33, 44, 46
P 6024	PUTRESCINE DYHROCHLORIC ACID	1.0 mg	To prepare 161.1 µg/ml (1mM) stock solution; add 6.2 ml sterile medium, gently swirl to dissolve.	16.1 µg/ml (100 mM)	powder-RT; stock-frozen in working aliquots, avoid repeated freeze/thaw; solution stable 2-8°C, 30 days	25, 48

**Defined Medium Supplements** continued

S 9133	SODIUM SELENITE	1.0 mg	To prepare 20 µg/ml stock solution; add 50 ml sterile culture medium.	≤0.2 µg/ml	powder-RT; stock-frozen in working aliquots, avoid repeated freeze/thaw; solution stable 2-8°C, 30 days	7, 24, 40, 41, 50, 52
S 0885	SOMATOSTATIN	0.02 mg	To prepare 2 µg/ml stock solution; add 10 ml sterile culture medium.	0.3-50 ng/ml	powder -0°C stock-frozen in working aliquots, avoid repeated freeze/thaw	20
T 5641	TESTOSTERONE	1.0 mg	To prepare 20 µg/ml stock solution; add 2.0 ml abs. ethanol, gently swirl to dissolve, add 48 ml sterile medium.	20-300 ng/ml	powder-RT; stock-frozen in working aliquots, avoid repeated freeze/thaw, solution stable 2-8°C, 30 days	3, 9, 21, 50
T 4393	THROMBIN	0.1 mg	To prepare a 10 µg/ml stock solution; add 10 ml sterile culture medium.		powder -0°C; stock-frozen in working aliquots, avoid repeated freeze/thaw, solution stable 2-8°C, 30 days	6, 26
T 9146	THYROTROPIN-RELEASING HORMONE	1.0 mg	To prepare 20 µg/ml stock solution; add 50 ml sterile culture medium.	0.3-10 ng/ml	powder -0°C; stock-frozen in working aliquots, avoid repeated freeze/thaw, solution stable 2-8°C, 30 days	22, 37
T 0397	L-THYROXINE	1.0 mg	To prepare 20 µg/ml stock solution; add 50 ml sterile culture medium.	5-50 ng/ml	powder -0°C; stock-frozen in working aliquots, avoid repeated freeze/thaw; solution stable 2-8°C, 30 days	23, 38, 51
T 5391	TRANSFERRIN HUMAN	10 mg	To prepare 500 µg/ml stock solution; add 20 ml sterile tissue culture medium, gently swirl to dissolve.	0.5-100 µg/ml	powder 2-8°C stock/frozen in working aliquots, avoid repeated freeze/thaw, solution stable 2-8°C for 30 days	15, 48

## Defined Medium Supplements continued

T 5516	3,3',5-TRIIODO-L- THYRONINE	1.0 mg	To prepare 20 µg/ml stock solution; add 1.0 ml 1.0 N NaOH, gently swirl to dissolve, add 49 ml sterile medium.	0.02-50 ng/ml	powder -0°C; stock-frozen in working aliquots, avoid repeated freeze/thaw, solu- tion stable 2-8°C, 30 days	9, 28, 45
--------	--------------------------------	--------	--	---------------	--	-----------

## REFERENCES

1. Armanini, D. et.al. (1987). *Jour. of Clin. Endo. and Met.* 65:101-104.
2. Armanini, D. et.al. (1987). *Jour. of Steroid Biochem.* 27:967-969.
3. Barnes, D. and Sato, G. (1980). *Analytical Biochem.* 102:255-270.
4. Barnes, D. and Sato, G. (1980). *Cell* 22:649.
5. Bell, N. (1970). *J. Clin. Invest.* 49:1368.
6. Biggs, R. (1977). *Human Blood Coagulation, Haemostasis and Thrombosis, Second Edition.* Blackwell Scientific Publications, Philadelphia, PA. p.772.
7. Bottenstein, J. and Sato, G. (1978). *Proc. Natl. Acad. Sci. USA* 78:514.
8. Bridson, W. and Kohler, P. (1970). *J. Clin. Endocrinol. Metab.* 30:538.
9. Colowick, S. and Kaplan, N. (1979). *Methods in Enzymology.* Vol 58:99-103.
10. Fischer, G. and Wiesner, R. (1983). *Hormonally Defined Media* Springer-Verlag Berlin, Heidelberg, Germany. pp. 235,349.
11. *ibid.*, (1983). pp. 6, 80, 147, 345.
12. *ibid.*, (1983). pp. 134, 216, 224, 230, 234.
13. *ibid.*, (1983). pp. 330, 335, 340, 352, 368, 435.
14. *ibid.*, (1983). pp. 46, 79, 120, 123, 287.
15. *ibid.*, (1983). pp. 6, 19-25.
16. *ibid.*, (1983). p. 123.
17. *ibid.*, (1983). pp. 234, 264, 274.
18. *ibid.*, (1983). pp. 120-123.
19. *ibid.*, (1983). pp. 19, 58, 107, 120, 436.
20. *ibid.*, (1983). p. 286.
21. *ibid.*, (1983). p. 447.
22. *ibid.*, (1983). pp. 134, 163, 252, 294.
23. *ibid.*, (1983). pp. 190, 210, 285, 407.
24. *ibid.*, (1983). pp. 134, 144, 180, 190, 216, 222, 230, 252.
25. *ibid.*, (1983). pp. 6, 216, 252, 330.
26. *ibid.*, (1983). pp. 88, 90.
27. Guibert, L. and Iscove, N. (1976) *Nature (London)* 263:594.
28. Hayashi, I. et al. (1978). *In Vitro* 14:23.
29. Hirsch, P. and Munson, P. (1969). *Physiol. Rev.* 49:548.
30. Iscove, N. and Melchers, F. (1978). *J. Exp. Med.* 147:923.
31. Jimenez de Asua, L. et.al. (1979). *Hormones in Cell Culture.* Cold Spring Harbor Conferences in Cell Proliferation. Sato, G. and Ross, R. (eds). Cold Spring Harbor Laboratory, Cold Spring Harbor, NY. Vol. 6:403-424. Jimenez de Asua, L. et.al. (1975). *Proc. Natl. Acad. Sci. USA.* 72: 2724-2728.
33. Jimenez de Asua, L. et.al. (1982). *Prostaglandins and Cancer: First International Conference.* Alan R. Liss, Inc., New York, NY. pp. 309-331.
34. Kahwanago, I. et.al. (1970). *Endocrinology* 86:1319.
35. Kohler, P. et.al. (1968). *Biochem. Biophys. Res. Commun.* 33:834.
36. Kruse, P. and Patterson, K. (1973). *Tissue Culture Methods and Applications.* Academic Press, New York, NY. pp. 564-575.
37. Kruse, P. and Patterson, K. (1971). *Tissue Culture Methods and Applications.* Academic Press, New York, NY. pp. 584-586.

38. Lauder, J. (1978). Brain Res. 142:25-39.
39. Mather, J. and Sato, G. (1979). Exp. Cell Res. 120:191-200.
40. McKeehan, W. et al. (1976). Proc. Natl. Acad. Sci. USA. 73:2023.
41. Murakami, H. and Masui, H. (1980). Proc. Natl. Acad. Sci. USA. 77:3464.
42. Notides, A. (1971). Endocrinology 89:229.
43. Otto, A. et.al. (1982). Proc. Natl. Acad. Sci. USA. 79:4992-4996.
44. Powles, T. et.al. (1982). Prostaglandins and Cancer:First International Conference. Alan R. Liss, Inc., New York, NY.
45. Samuels, H. et al. (1973). Science 181:1253.
46. Samuelson, B. and Poletti, R. (1976). Advances in Prostaglandin and Thromboxane Research. Raven Press, New York, NY.
47. Sato, G. and Ross R. (1979). Hormones and Cell Culture, Book B. Cold Spring Harbor Laboratory. pp. 655-668.
48. *ibid.*, (1979). Book B, pp. 524-555.
49. *ibid.*, (1979). Book B, pp. 536, 677, 683, 937-945.
50. Sato, G. and Reid, L. (1978). International Rev. of Biochem., Biochem and Mode of Action of Hormones II, (Univ. Park Press, Baltimore) Vol. 20, pp 219-251
51. Seelig, S. et al. (1981). Proc. Natl. Acad. Sci. USA. 78:4733-4737.
52. Stadtman, T. (1974). Science 183:915.
53. Waymouth, C. (1984). Preparation and Use of Serum-Free Culture Media. In: Methods for Preparation of Media, Supplements and Substrata For Serum-Free Animal Cell Culture. Alan R. Liss, Inc., New York, NY. Vol.1:23.
54. Wehling, M. et.al. (1987). ACTA Endocrinologica 116:555-560.
55. Wehling, M. et.al. (1987). Amer. Jour. of Phys. 252:505-508.
56. Wehling, M. et.al. (1988). Clin. Endo. (Oxford). 28:67-74.