


# MultiScreen<sup>®</sup> Filter Plate with Ultracel<sup>®</sup>-10 Membrane

protocol note

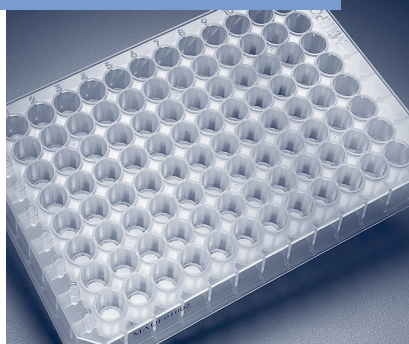

---

## ***Protein Retention, Recovery, Volume Recovery, and Guidelines for Concentration and Desalting***

### **Introduction**

MultiScreen Filter Plate with Ultracel-10 membrane (MultiScreen Ultracel-10 filter plate) provides a new method for high throughput sample preparation. The ultrafiltration-based filter plate is designed for automation-compatible sample purification, concentration and desalting of biological solutions, as well as protein removal from samples prior to instrument analysis. The 96-well MultiScreen filter plate incorporates the Ultracel-10 membrane (10,000 nominal molecular weight limit regenerated cellulose) for ultra low-binding, high recovery results. It is designed for use with centrifugation.

This protocol note includes guidelines for protein retention, recovery and concentration/desalting using MultiScreen Ultracel-10 filter plates in centrifugal mode. Ultrafiltration is shown to be a highly effective and rapid means for these applications. The 96-well filter plate increases sample throughput and is compatible with standard microtiter plates, instrumentation and liquid handling equipment.

Part 1: **Protein Retention**

Part 2: **Protein Recovery**

Part 3: **Protein Desalting**

Part 4: **Ultrafiltration Time and Volumes**

## Part 1: Protein Retention

**Objective:** The goal of this experiment was to determine the protein retention performance, as well as ultrafiltration reliability across the plate as measured by protein recovery. FITC-BSA (67 kDa) and Cytochrome c (12 kDa) were used as model molecules in this experiment.

### Materials

#### Reagents

- Milli-Q® water
- Phosphate Buffered Saline (PBS) (Sigma-P 3813, pH 7.4)
- Cytochrome c (Sigma C-8266), 0.5 mg/mL solution in PBS
- Bovine Albumin – Fluorescein Isothiocyanate labeled (FITC-BSA) (SIGMA A-9771), 0.5 mg/mL solution in PBS

#### Equipment and Materials

- Centrifuge equipped with a 96-well plate rotor
- Molecular Devices 96-well plate reader set to 410 nm wavelength. Reader must be equipped with auto mixer and Pathcheck™.
- SOFTMax® Pro 3.1 software for Molecular Devices reader (Reader must be equipped with auto mixer and Pathcheck).
- Wallac/Victor™ Fluorescence Detector or comparable instrument
- 96-well collection plates (Costar 9017 or equivalent)
- Black 96-well collection plates (Costar 3915 or equivalent)
- Plate shaker


### Protocol (Part 1 Protein Retention)

1. Place a MultiScreen Ultracel-10 filter plate on top of a 96-well collection plate (Costar 9017).
2. Add 200  $\mu\text{L}$ /well of Cytochrome c or 300  $\mu\text{L}$ /well of FITC-BSA to the MultiScreen Ultracel-10 filter plate.
3. Centrifuge at 3000 x g for 1 hour, until wells are empty.
4. Evaluate the concentration of protein in the filtrate as follows.
  - a. Determine FITC-BSA concentration by reading plates at 485Ex and 535Em using a Fluorescence Detector.
  - b. Determine Cytochrome c concentration by reading plates at 410 nm wavelength using a spectrophotometer.
5. After generating a standard curve for both proteins, calculate protein concentrations in both the retentates and filtrates from the standard curve. Calculate percent protein retention based on the amount of protein in the filtrate divided by the initial protein concentration of 0.5mg/mL.


### Results

The results show average Cytochrome c retention greater than 97% and average BSA retention greater than 99.9% (limit) with excellent well-to-well reproducibility (see Figures 1 and 2).

**Figure 1.** Average protein retention of Cytochrome c (n = 76 per well position, sd <10%).

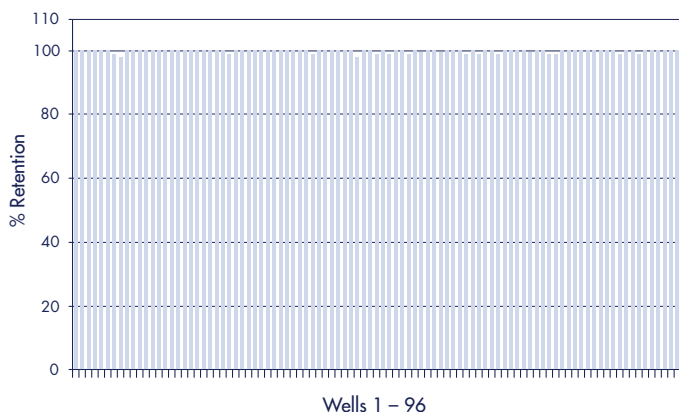

**Figure 2.** Average protein retention for BSA-FITC (n = 76 per well position, sd <10%).

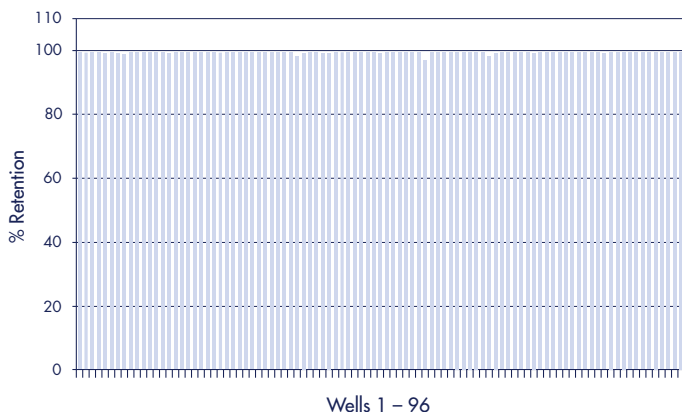

## Part 2: Protein Recovery

**Objective:** Ultrafiltration provides an effective means for the rapid concentration of biological fluids and protein solutions. With Ultracel-10 ultrafiltration membrane, it is expected that any protein over 10 kDa will be retained by the membrane. The goal of this experiment was to determine protein recovery, as well as ultrafiltration reproducibility across the plate as measured by protein recovery. FITC-BSA (67 kDa) and Cytochrome c (12 kDa) were used as a model molecules in this experiment.

### Method

#### Reagents

- Milli-Q water
- Phosphate Buffered Saline (PBS) (Sigma P 3813, pH 7.4)
- Cytochrome c (Sigma C-8266), 0.5 mg/mL solution in PBS
- Bovine Albumin – Fluorescein Isothiocyanate labeled (FITC-BSA) (SIGMA A-9771), 0.5 mg/mL solution in PBS

#### Equipment and Materials

- Centrifuge equipped with 96 well plate rotor
- Molecular Devices 96 well plate reader set to 410 nm wavelength. Reader must be equipped with auto mixer and Pathcheck.
- SOFTMax Pro 3.1 software for Molecular Devices reader
- Wallac/Victor Fluorescence Detector or comparable instrument
- 96-well collection plates (Costar 9017 or equivalent)
- Black 96-well collection plates (Costar 3915 or equivalent)
- Plate shaker

#### Protocol (Part 2 Protein Recovery)

1. Place the MultiScreen Ultracel-10 filter plate on top of the 96-well collection plate (Costar 9017).
2. Add 200  $\mu\text{L}$ /well of Cytochrome c or 300  $\mu\text{L}$ /well of FITC-BSA to the MultiScreen Ultracel-10 filter plate.
3. Centrifuge at 3000  $\times$  g for 1 hour, until wells are empty.
4. Add 200  $\mu\text{L}$ /well or 300  $\mu\text{L}$ /well PBS to each well with a multi-channel pipettor.
5. Agitate the plates for 3 minutes on a plate shaker.
6. Transfer the resuspended retentate into a clear 96-well collection plate for Cytochrome c and a black collection plate for FITC-BSA analysis.
  - a. Determine FITC-BSA concentration by reading plates at 485Ex and 535Em using a Fluorescence Detector.
  - b. Determine Cytochrome c concentration by reading plates at 410 nm wavelength using a spectrophotometer.
7. After generating a standard curve for both proteins calculate protein concentrations in both the retentate and filtrates from the standard curve. Calculate percent protein recovery based on the initial protein concentration of 0.5 mg/mL.

**Table 1.** Average protein recovery for Cytochrome c and FITC-BSA after concentration in MultiScreen Ultracel-10 filter plates (n=96).


	<b>Cytochrome c</b>	<b>FITC-BSA</b>
Plate 1	85%	88%
Plate 2	92%	87%

## Results


Table 1 shows the results of protein recovery for 200  $\mu\text{L}$  of 0.5 mg/mL of Cytochrome c and FITC-BSA, based on 2 plates (192 wells) analyzed for each protein. The protein solutions were spun to dryness and resuspended in 50  $\mu\text{L}$  of PBS. Figure 3 shows individual recovery numbers for 96 wells tested with Cytochrome c. The results demonstrate that over 85% of either protein can be recovered after 1 hour spin-to-dryness, and that recovery is consistent throughout all 96 wells of MultiScreen Ultracel-10 filter plates.

Even higher recoveries of the protein can be achieved with increased resuspension volume and time on a shaker. Figure 4 shows that 98.5% of protein can be recovered if the protein is resuspended in 200  $\mu\text{L}$ , and the plate is agitated for 30 minutes for resuspension.

**Figure 3.** Percent retentate starting from 0.5 mg/mL Cytochrome c.

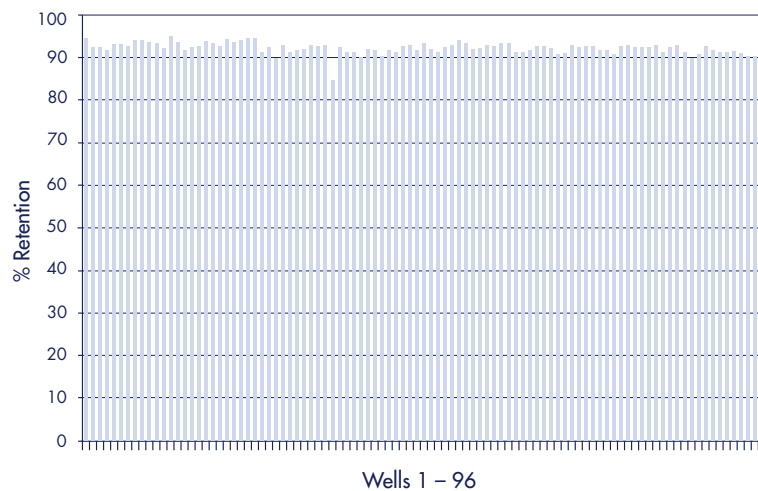

**Figure 4.** The effect of resuspension volume and shake time.

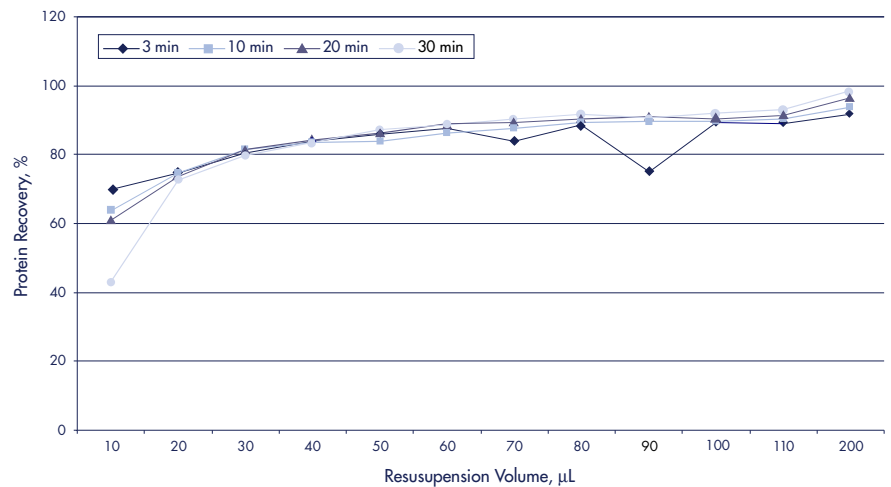

**Table 2.** Well-to-well reproducibility of Cytochrome c recovery.

avg % retentate	92.5	min	84.9
std dev	1.3	max	95.4

## Part 3: Protein Desalting

**Objective:** The MultiScreen Ultracel-10 filter plates is an efficient tool for the removal of salts, unincorporated radioisotopes, dye and other small molecules from protein solutions. Usually it requires more than one round of ultrafiltration to remove a majority of salt and exchange protein into low salt buffer. To achieve this, concentrated protein gets re-dissolved in water or low molarity buffer and concentrated a second time. Multiple dilution/concentration steps may be required until salt is removed. It is important that proteins be recovered from the membrane after repeated centrifugations. In this experiment, we demonstrate MultiScreen Ultracel-10 filter plate performance for desalting of proteins.

### Method

#### Reagents

- Milli-Q water
- Phosphate Buffered Saline (PBS) (Sigma P 3813, pH 7.4)
- Cytochrome c (Sigma C-8266), 0.5 mg/mL solution in PBS
- Bovine Albumin – Fluorescein Isothiocyanate labeled (FITC-BSA) (SIGMA A-9771), 0.5 mg/mL solution in PBS

#### Equipment and Materials

- Centrifuge equipped with 96-well plate rotor
- Molecular Devices 96 well plate reader set to 410 nm wavelength. Reader must be equipped with auto mixer and Pathcheck.
- SOFTMax Pro 3.1 software for Molecular Devices reader
- 96-well collection plates (Costar 9017 or equivalent)
- Black 96-well collection plates (Costar 3915 or equivalent)
- Plate shaker
- Conductivity meter


#### Protocol (Part 3 Protein Desalting)

1. Prepare 0.25 mg/mL solution of Cytochrome c in 0.5M NaCl.
2. Assemble three MultiScreen Ultracel-10 filter plates on top of 96-well collection plates.
3. Add 200  $\mu$ L/well of protein solution to the MultiScreen Ultracel-10 filter plate.
4. Centrifuge at 3000 x g for 1 hour, until wells are empty
5. After the MultiScreen Ultracel-10 filter plate has spun, add 200  $\mu$ L water to each well.
6. Shake the MultiScreen Ultracel-10 filter plate for 3 minutes.
7. Pool the protein solution from 96 wells of one plate together. Measure conductivity and determine protein concentration at 410 nm wavelength.
8. Repeat steps 4 – 7 for remaining two plates. Analyze the second plate after the second cycle and analyze the third plate after the third cycle.

**Table 3.** Salt removal and protein recovery using the MultiScreen Ultracel-10 filter plate. The starting solution was 0.25 mg/mL of Cytochrome c in 0.5M NaCl.

	% Protein Recovery	% Salt Remaining
Spin 1	92.8%	5%
Spin 2	89.7%	0.3%
Spin 3	85.8%	0.04%

**Figure 5.** Percentage of Cytochrome c and NaCl remaining in the Ultracel-10 wells after one, two, or three concentration/dilution cycles.

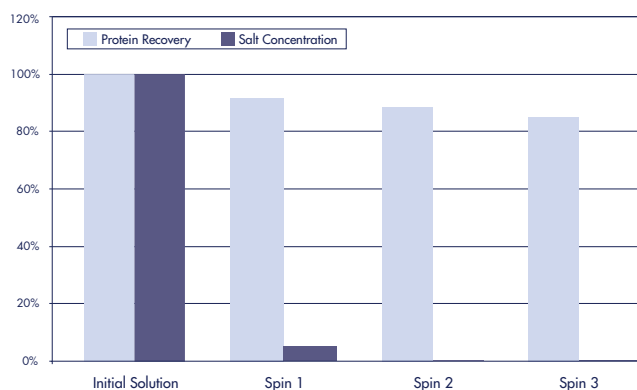

### Results

Table 3 and Figure 5 show the results of salt removal and protein recovery in MultiScreen Ultracel-10 filter plates after 3 consecutive centrifugations of Cytochrome c solution with 0.5M NaCl. It is demonstrated that up to 95% of salt can be removed in the first

cycle of concentration/desalting. As much as 99.9% salt can be removed after two centrifugations. It is important to note that protein recovery numbers are still high, with around 85% of protein being retained by the membrane after three centrifugation cycles.

## Part 4: Ultrafiltration Time and Volumes

**Objective:** In this experiment, the time required to spin samples to dryness as a function of starting volume and protein concentration was determined.

### Method

#### Reagents

- Milli-Q water
- Phosphate Buffered Saline (PBS) (Sigma P 3813, pH 7.4)
- Cytochrome c (Sigma C-8266), 0.5 mg/mL solution in PBS
- Bovine Albumin – Fluorescein Isothiocyanate labeled (FITC-BSA) (SIGMA A-9771), 0.5 mg/mL solution in PBS


#### Equipment and Materials

- Centrifuge equipped with a 96-well plate rotor
- Molecular Devices 96-well plate reader set to 410 nm wavelength. Reader must be equipped with auto mixer and Pathcheck.
- SOFTMax Pro 3.1 software for Molecular Devices reader
- Wallac/Victor Fluorescence Detector or comparable instrument
- 96-well collection plates (Costar 9017 or equivalent)
- Black 96-well collection plates (Costar 3915 or equivalent)
- Plate shaker

#### Protocol

1. Prepare 0.5 mg/mL, 5 mg/mL and 10 mg/mL solutions of Cytochrome c in PBS.
2. Place a MultiScreen Ultracel-10 filter plate on top of a 96-well collection plate.
3. Add 100  $\mu\text{L}$ /well, 200  $\mu\text{L}$ /well and 300  $\mu\text{L}$ /well of protein solution to different wells of a MultiScreen Ultracel-10 filter plate.
4. Place plates in centrifuge and spin at 3000 x g. Check for retentate volumes at 15, 25, 30, 45, 60, 75, 90, 105 and 150 minutes.

**Figure 6.** Times required to spin 3 concentrations of Cytochrome c solutions to dryness at 3000 x g, 25 °C.

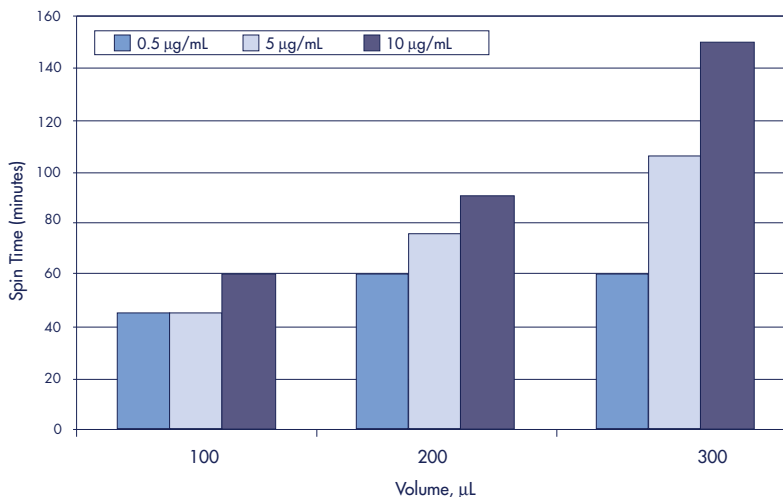


## Results

Results of this study are shown in Table 4 and Figure 6. The time to dryness depends on the protein concentration and starting volume. As long as 150 minutes centrifugation may be required to concentrate 300  $\mu\text{L}$  of 10 mg/mL Cytochrome c solution. Figure 7 demonstrates that volume reduction is reproducible throughout all 8 rows of wells of the MultiScreen Ultracel-10 filter plate.

**Table 4.** Times required to spin to 50  $\mu\text{L}$  retentate at 3000 x g, 25 °C

Initial Volume of 0.5 mg/mL Cytochrome c ( $\mu\text{L}$ )	Time to Spin to 50 $\mu\text{L}$ Retentate (minutes)
100 $\mu\text{L}$	15
200 $\mu\text{L}$	25
300 $\mu\text{L}$	30

**Figure 7.** Row-to-row variations of retentate volumes for Cytochrome c at various times at 25 °C and 3000 x g with 300  $\mu\text{L}$  starting volume.

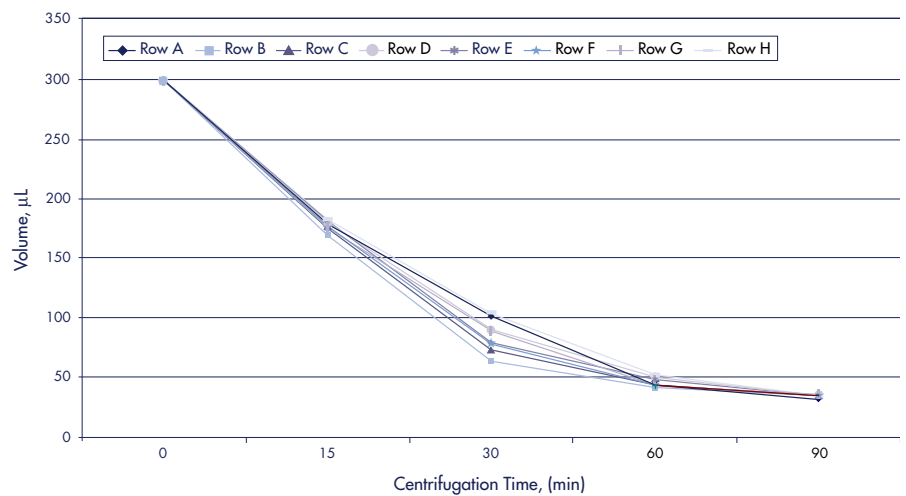

## Summary and Conclusions

Protein retention, recovery, volume recovery, concentration and desalting can be accomplished rapidly and with good results using centrifugation and an ultrafiltration membrane. Results presented here and published separately demonstrate that MultiScreen Filter Plate with Ultracel-10 membrane is designed for high retention and low protein binding. It is an efficient tool for the parallel processing of 96 samples at once by centrifugation.

## Related literature

PF2050EN00: MultiScreen Filter Plate with Ultracel-10 Membrane data sheet

## To Place an Order or Receive Technical Assistance

For additional information call your nearest Millipore office:  
In the U.S. and Canada, call toll-free **1-800-MILLIPORE (1-800-645-5476)**  
In the U.S., Canada and Puerto Rico, fax orders to **1-800-MILLIFX (1-800-645-5439)**  
Internet: [www.millipore.com](http://www.millipore.com)  
Technical Service:  
[www.millipore.com/techservice](http://www.millipore.com/techservice)



Now you can buy Millipore products online @ [www.millipore.com/purecommerce](http://www.millipore.com/purecommerce)

Millipore, MultiScreen, Ultracel and Milli-Q are registered trademarks of Millipore Corporation.  
Pathcheck and SoftMax are registered trademarks of Molecular Devices CCorproation.  
Wallac Victor is a trademark of PerkinElmer or an affiliated company.  
Lit No. PC1025EN00 8/03 03-235  
© 2003 Millipore Corporation, Billerica, MA 01821  
U.S.A. All rights reserved.

## Millipore Worldwide

### AUSTRALIA

Tel. 1 800 222 111  
or (02) 9888 8999  
Fax (02) 9878 0788

### AUSTRIA

Tel. 0820 87 44 64  
Fax 0820 87 44 44

### BALTIC COUNTRIES

Tel. +358 203 05 645  
Fax +358 203 05 644

### BELGIUM

Tel. 070 225 645  
Fax 070 225 644

### BRAZIL

Tel. (011) 5548-7011  
Fax (011) 5548-7923

### CANADA

Tel. (800) 645-5476  
Fax (800) 645-5439

### CHINA, PEOPLE'S REPUBLIC OF

Beijing:  
Tel. (86-10) 8519 1250  
(86-10) 6518 1058  
Fax (86-10) 8519 1255  
Guangzhou:  
Tel. (86-20) 8752 0187  
(86-20) 8752 0173  
Fax (86-20) 8752 0172  
Hong Kong:  
Tel. (852) 2803 9111  
Fax (852) 2513 0313  
Shanghai:  
Tel. (86-21) 5306 9100  
Fax (86-21) 5306 0838

### CZECH REPUBLIC

Tel. 02 2051 3841  
Fax 02 2051 4298

### DENMARK

Tel. 7010 5645  
Fax 7010 5644

### EASTERN EUROPE, C.I.S., AFRICA, MIDDLE EAST AND GULF

Tel. +33 3 88 38 9536  
Fax +33 3 88 38 9539

### FINLAND

Tel. 0203 05 645  
Fax 0203 05 644

### FRANCE

Tel. 0825 045 645  
Fax 0825 045 644

### GERMANY

Tel. 01 805 045 645  
Fax 01 805 045 644

### HUNGARY

Tel. 01 209 04 33  
Fax 01 209 02 95

### INDIA

Tel. 080 839 46 57  
Fax 080 839 63 45

### IRELAND

Tel. +44 1923 816375  
Fax +44 1923 818297

### ITALY

Tel. 848 845 645  
Fax 848 845 644

### JAPAN

Tel. (03) 5442-9711  
Fax (03) 5442-9736

### KOREA

Tel. (822) 551-0230  
Fax (822) 551-0228

### LUXEMBOURG

Tel. +32 2726 88 40  
Fax +32 2726 98 84

### MALAYSIA

Tel. 03-7957-1322  
Fax 03-7957-1711

### MEXICO

Tel. (55) 5576 9688  
Fax (55) 5576 8706

### THE NETHERLANDS

Tel. 0900 7 645 645  
Fax 0900 7 645 644

### NORWAY

Tel. 810 62 645  
Fax 810 62 644

### POLAND

Tel. 22 669 12 25  
Fax 22 663 70 33

### PORTUGAL

Tel. +34 91 728 39 60  
Fax +34 91 729 29 09

### PUERTO RICO

Tel. (787) 273-8495  
Fax (787) 747-6553

### SINGAPORE

Tel. 6842-1822  
Fax 6842-4988

### SPAIN

Tel. 901 516 645  
Fax 902 011 644

### SWEDEN

Tel. 0771 200 645  
Fax 0771 200 644

### SWITZERLAND

Tel. 0848 645 645  
Fax 0848 645 644

### TAIWAN

Tel. 886-2-2792-9333  
Fax 886-2-2792-6555

### U.K.

Tel. 0870 900 46 45  
Fax 0870 900 46 44

### U.S.A.

Tel. (800) 645-5476  
Fax (800) 645-5439

### IN ALL OTHER COUNTRIES

Millipore International  
Tel. +1 781 533 8622  
Fax +1 781 533 8630

# MILLIPORE