

Product Information

EPIDERMAL GROWTH FACTOR (EGF)

Tissue Culture Grade
From Mouse Submaxillary Glands

Product Number **E4127**

Product Description

Epidermal Growth Factor (EGF) is a 6 kDa polypeptide discovered by Cohen and Levi-Montalcini and isolated from mouse submaxillary glands.^{1,2,3} EGF is mitogenic for a variety of epidermal and epithelial cells, including fibroblasts, glial cells, mammary epithelial cells, vascular and corneal endothelial cells, bovine granulosa, rabbit chondrocytes, HeLa and SV40-3T3 cells.⁴ In tissue cultures EGF can act to reduce or eliminate the requirement for serum and is often used in conjunction with other media additives and hormones, such as Insulin (Product No. I1882), Transferrin (Product No. T5391), and Prostaglandin E₁ (Product No. P7527).⁵ Cellular metabolic effects of EGF include stimulation of ion fluxes,⁶ glucose transport,⁷ glycolysis,⁸ and synthesis of DNA, RNA and proteins.⁴

Human EGF was originally isolated from human urine as beta-Urogastrone (Product No. U7126), an inhibitor of gastric acid secretion.⁹ Mouse and human EGF are 70% homologous in their sequence of 53 amino acids, act upon the same EGF receptor, elicit nearly identical biological effects and are highly inter-species cross-reactive.^{10,11,12,13}

EGF is structurally homologous to human Transforming Growth Factor-Alpha (Product Nos. T7533 and T2035), which exerts its actions through EGF receptors.¹⁴ EGF is also homologous to a sequence contained in a 19 kDa protein of Vaccinia virus,¹⁵ which appears to utilize the EGF receptor to gain entry into cells.¹⁶ The EGF receptor is a 170 kDa glycoprotein having EGF-activated protein tyrosine kinase activity.¹⁷ Platelet-Derived Growth Factor (PDGF, Product No. P8147) transmodulates the EGF receptor by reducing both its EGF affinity and its kinase activity.¹⁸

Human EGF is found in many body fluids, including urine, milk, saliva, sweat and seminal fluid,^{19,20} and has its highest concentration in alpha-granules of blood platelets.²¹ The biological role of EGF includes

the inhibition of gastric acid secretion,¹⁰ support of growth and differentiation during fetal development,²² neuromodulation in the central nervous system,²³ and stimulation of epidermal growth and keratinization.³ Clinical trials have shown that topical administration of human recombinant EGF accelerates wound healing.²⁴

Reagents

Lyophilized from a solution in 1 ml of 5 mM ammonium acetate, pH 6.5.

Storage/Stability

Store vial at 2-8 °C.

After reconstitution, the product may be stored for two weeks at 2-8 °C or may be stored as aliquots at -20 °C. Prolonged storage and/or repeated freezing and thawing of reconstituted product are not recommended.

Product Profile

To prepare a stock solution, reconstitute contents of the vial in a solution that contains 0.1-1.0% BSA or 1-10% serum in buffered saline or tissue culture media. This may be diluted before use to the final working concentration of EGF, generally 0.1-10 ng/ml. Additional filtration of the stock solution is not recommended and may result in product loss due to adsorption onto the filter membrane.

References

1. Cohen, S., Proc. Natl. Acad. Sci. USA, **46**, 302 (1960).
2. Levi-Montalcini, R., and Cohen, S., Ann. NY Acad. Sci, **85**, 324 (1960).
3. Cohen, S., J. Biol. Chem., **237**, 1555 (1962).
4. Carpenter, G. and Cohen, S., Annu. Rev. Biochem., **48**, 193 (1979).
5. Fischer, G. and Wieser, R., Hormonally Defined Media (Springer-Verlag Berlin, Heidelberg, Germany), p. 23 (1983).
6. Rozengurt, E. and Heppel, L. A., Proc. Natl. Acad. Sci. USA, **72**, 4492 (1975).

7. Barnes, D., and Colowick, S., J. Cell. Physiol., **89**, 633 (1976).
8. Diamond, I., et al., J. Biol. Chem., **253**, 866 (1978).
9. Cohen, S., and Carpenter, G., Proc. Natl. Acad. Sci. USA, **72**, 1317 (1975).
10. Gregory, H., Nature, **257**, 325 (1975).
11. Hollenberg, M., and Gregory, H., Life Sci., **20**, 267 (1977).
12. Carpenter, G., and Zendegui, J. G., Exp. Cell Res., **164**, 1 (1986).
13. George-Nascimento, C., et al. Biochemistry, **27**, 797 (1988).
14. Todaro, G. J., et al., Proc. Natl. Acad. Sci. USA, **77**, 5258 (1980).
15. Blomquist, M. C., et al., Proc. Natl. Acad. Sci. USA, **81**, 7363 (1984).
16. Eppstein, D. A., et al., Nature, **318**, 663 (1985).
17. Schlessinger, J., Biochemistry, **27**, 3119 (1988).
18. Davis, R. and Czech, M., J. Biol. Chem., **262**, 6832 (1988).
19. Hirata, Y. and Orth, D. N., J. Clin. Endocrinol. Metab., **48**, 673 (1979).
20. Oka, Y. and Orth, D. N., J. Clin. Invest. **72**, 249 (1983).
21. Carpenter, G. L., J. Cell Sci., Suppl., **3**, 1 (1985).
22. Thorburn, G., et al., Growth and Maturation Factors, Vol. 3 (G. Guroff, ed.) Wiley and Sons, NY, p. 175 (1986).
23. Fallon, J., et al., Science, **224**, 1107 (1984).
24. Brown, G., et al., N. Engl. J. Med., **321**, 76 (1989).
25. Carpenter, G. and Zendegui, J., Anal. Biochem., **153**, 279 (1985).
26. Taylor, J. M., et al., J. Biol. Chem., **247**, 5928 (1972).

JWM/DAA 2/2003

Sigma brand products are sold through Sigma-Aldrich, Inc.

Sigma-Aldrich, Inc. warrants that its products conform to the information contained in this and other Sigma-Aldrich publications. Purchaser must determine the suitability of the product(s) for their particular use. Additional terms and conditions may apply. Please see reverse side of the invoice or packing slip.