

Product Information

2,2'-Azino-bis(3-ethylbenzothiazoline-6-sulfonic acid) diammonium salt tablet

10 mg substrate per tablet

A9941

Product Description

CAS Number: 30931-67-0 [2,2'-Azino-bis(3-ethylbenzothiazoline-6-sulfonic acid) diammonium salt component]


Molecular Formula: $C_{18}H_{24}N_6O_6S_4$ [2,2'-Azino-bis(3-ethylbenzothiazoline-6-sulfonic acid) diammonium salt component]

Molecular Weight: 548.68 [2,2'-Azino-bis(3-ethylbenzothiazoline-6-sulfonic acid) diammonium salt component]

Synonym: AzBTS-(NH_4)₂, ABTS™-(NH_4)₂, Diammonium 2,2'-azino-bis(3-ethylbenzothiazoline-6-sulfonate)

Storage Temperature: Room temperature

Structure [2,2'-Azino-bis(3-ethylbenzothiazoline-6-sulfonic acid) diammonium salt component]:

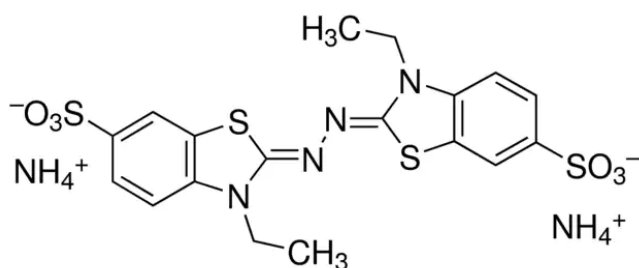

2,2'-Azino-bis(3-Ethylbenzthiazoline-6-Sulfonic Acid), also known as AzBTS, is a chromogen that is suitable for use in ELISA procedures which utilize horseradish peroxidase (HRP) conjugates.¹⁻³ The reaction of the ABTS substrate with HRP produces a soluble end product that is green in color and can be read spectrophotometrically at 405 nm.⁴ Several theses^{5,6} and dissertations⁷⁻¹³ cite use of this A9941 tablet product in their protocols.

Precautions and Disclaimer

For R&D use only. Not for drug, household, or other uses. Please consult the Safety Data Sheet for information regarding hazards and safe handling practices.

Reagent

Tablet Properties

- Substrate Content: 10 mg
 - Appearance: Green, round tablet, 5/32 in. diameter
 - Tablet Weight: 35 mg (range 31.5-38.5 mg)
 - Solubility (*): Clear, colorless to a clear, light-green solution
 - Dissolution Time (*): Not more than 5.0 minutes
 - pH: 5.0 (range 4.2-5.8) in 100 mL of deionized water
 - Activity: 100% (range 90-110%)
 - Packaging: 50 or 100 tablets per box, individually foil wrapped for ease of use, storage, and safety
- (*): One tablet is dissolved in 100 mL of deionized water.

Preparation Instructions

Option 1

- Dissolve 1 tablet in 100 mL of 0.05 M phosphate-citrate buffer, pH 5.0 (such as Cat. No. P4809).
- Add 25 μ L of fresh 30% hydrogen peroxide (H_2O_2 , such as Cat. No. H1009) immediately prior to use.

Option 2

- Dissolve 1 tablet in 100 mL of 0.05 M phosphate-citrate buffer, pH 5.0, containing 0.03% sodium perborate (such as Cat. No. P4922).

Phosphate-Citrate Buffer Preparation

To prepare 0.05 M phosphate-citrate buffer, pH 5.0:

1. Add 25.7 mL of 0.2 M dibasic sodium phosphate (such as Cat. Nos. S0876 or 71643) to 24.3 mL of 0.1 M citric acid (such as Cat. Nos. C7129 or C0706).
2. Adjust the buffer volume to 100 mL with deionized water.
3. Adjust the pH to 5.0, if necessary.

References

1. Jones, A.R. *et al.*, *Biotechnol. J.*, **9(5)**, 664-674 (2014).
2. Yadirgi, G. *et al.*, *J. Immunol. Methods*, **451**, 90-99 (2017).
3. Gallati, V.H., *J. Clin. Chem. Clin. Biochem.*, **17(1)**, 1-7 (1979).
4. Porstmann, B. *et al.*, *J. Clin. Chem. Clin. Biochem.*, **19(7)**, 435-439 (1981).
5. Custers, Jean, "Identifying the gene involved in the shape of potato tubers". Wageningen University, M.Sc. thesis, p. 18 (2015).
6. Tardón, Daniela Irlanda Castro, "Integration between healthy and microbiome in Holstein cows immediately after calving". Universidade de São Paulo, M.Sc. thesis, p. 47 (2020).
7. Driedger, Darcy Renard, "Analysis of Potato Glycoalkaloids by Immunoassay Coupled to Capillary Electrophoresis or Matrix-Assisted Laser Desorption/Ionization Mass Spectrometry". University of Alberta, Ph.D. dissertation, pp. 106, 117 (2000).
8. Seymour, Maile Elizabeth Kahawaluaakalani, "Process Considerations in the Extraction and Recovery of Plant Virus Vaccines". University College London, Ph.D. dissertation, p. 38 (2001).
9. Li, Keyu, "Streamlined Development of Fully Human Antibody Fragments for ImmunoPET Imaging Using Phage Display Technology". University of California Los Angeles, Ph.D. dissertation, p. 38 (2014).

10. McGiven, John, "Immunoselection and structural evaluation of *Brucella* O-polysaccharide epitopes and their application to the serodiagnosis of bovine brucellosis". Imperial College London, Ph.D. dissertation, p. 140 (2014).
11. Lehmann, Sabrina, "Biology of Odoriferous Defensive Stink Glands of the Red Flour Beetle *Tribolium castaneum*". Georg-August-Universität Göttingen, Dr. rer. nat. dissertation, p. 27 (2015).
12. Otieno, Brunah Akinyi, "Semi-automated Ultrasensitive Electrochemical Microfluidic Immunoarray for On-line Capture and Detection of Cancer Protein Biomarkers". University of Connecticut, Ph.D. dissertation, p. 145 (2016).
13. Frentzel, Sarah, "IkB_{NS} differentially affects *in vivo* CD4⁺ and CD8⁺ T cell activation and plays a detrimental role in innate immunity to *Listeria monocytogenes* infection in mice". Otto-von-Guericke-Universität Magdeburg, Dr. rer. nat. dissertation, p. 21 (2018).

Notice

We provide information and advice to our customers on application technologies and regulatory matters to the best of our knowledge and ability, but without obligation or liability. Existing laws and regulations are to be observed in all cases by our customers. This also applies in respect to any rights of third parties. Our information and advice do not relieve our customers of their own responsibility for checking the suitability of our products for the envisaged purpose.

The information in this document is subject to change without notice and should not be construed as a commitment by the manufacturing or selling entity, or an affiliate. We assume no responsibility for any errors that may appear in this document.

Technical Assistance

Visit the tech service page at SigmaAldrich.com/techservice.

Terms and Conditions of Sale

Warranty, use restrictions, and other conditions of sale may be found at SigmaAldrich.com/terms.

Contact Information

For the location of the office nearest you, go to SigmaAldrich.com/offices.

The life science business of Merck KGaA, Darmstadt, Germany operates as MilliporeSigma in the U.S. and Canada.

MilliporeSigma, and Sigma-Aldrich are trademarks of Merck KGaA, Darmstadt, Germany or its affiliates. All other trademarks are the property of their respective owners. Detailed information on trademarks is available via publicly accessible resources.

© 1999-2022 Merck KGaA, Darmstadt, Germany and/or its affiliates. All Rights Reserved.

A9941dat Rev 10/22 PCS,GCY,MAM

**MILLIPORE
SIGMA**