



3050 Spruce Street
Saint Louis, Missouri 63103 USA
Telephone 800-325-5832 • (314) 771-5765
Fax (314) 286-7828
email: techserv@sial.com
sigma-aldrich.com

Product Information

Phenazine methosulfate

Product Number **P 9625**
Storage Temperature -0 °C

Product Description

Molecular Formula: $C_{14}H_{14}N_2O_4S$
Molecular Weight: 306.3
CAS Number: 299-11-6
Extinction coefficient: E^M for oxidized phenazine methosulfate (PMS) at maximum wavelength 387 nm is 26,300 (pH 2-8). The difference in molar extinction coefficients between oxidized and reduced PMS is 25,000.¹
Melting Point: 158-160 °C

PMS is an electron acceptor and carrier in enzyme systems. The oxidized form is yellow and the reduced form is colorless. Since the reduced PMS is easily oxidized by oxygen, it is used in assays as an electron carrier between enzymes and oxygen, cytochrome c, indophenols, or tetrazolium salts. The reduced PMS is used as an electron donor to reduce cytochrome c or in photosynthetic experiments. PMS is reduced by flavoproteins such as succinic dehydrogenase. PMS is reduced non-enzymatically by NADH and NADPH. PMS is also reduced by dithionite, sodium borohydride, ascorbic acid, reduced ubiquinones, and reduced vitamin K. PMS reacts with sulfhydryl groups including those of enzymes.¹ Usage of PMS for detection of specific dehydrogenases has been reported.^{2,3}

Precautions and Disclaimer

For Laboratory Use Only. Not for drug, household or other uses.

Preparation Instructions

Phenazine methosulfate is soluble in water (approximately 200 mg/ml).

Storage/Stability

Solutions should be prepared in deionized water, and not in neutral buffers. Solutions in water are stable frozen and protected from light for several months. Decomposition of solutions (with formation of pyocyanine and other products) may take hours when they are exposed to diffuse light, but only 5-10 minutes when exposed to sunlight.

Solutions are stabilized by polyvalent cations. Cations such as Ca^{2+} , Mg^{2+} , Mn^{2+} , Co^{2+} , Cu^{2+} , Ba^{2+} , Zn^{2+} , and Sn^{3+} at 100 mM each, prevented decomposition of PMS solutions when stored for 1 month, but were not protective when solutions were exposed to sunlight.^{1,2}

References

1. Data for Biochemical Research, 3rd ed., Dawson, R. M. C., et al., Oxford University Press (New York, NY: 1986), pp. 356-357.
2. Nachlas, et al., J. Biol. Chem., **235**, 499-503 (1960).
3. Babson, A. L., and Babson, S. R., Clin. Chem., **19(7)**, 766-769 (1973).

ARO/RXR 9/02

Sigma brand products are sold through Sigma-Aldrich, Inc.

Sigma-Aldrich, Inc. warrants that its products conform to the information contained in this and other Sigma-Aldrich publications. Purchaser must determine the suitability of the product(s) for their particular use. Additional terms and conditions may apply. Please see reverse side of the invoice or packing slip.

