

Product Information

Apyrase from potato

Recombinant, expressed in *Pichia pastoris*, ATPase ≥ 1000 units/mg protein, lyophilized powder

A6237


Product Description

CAS Number: 9000-95-7

Enzyme Commission (EC) Number: 3.6.1.5

Synonyms: Nucleoside-triphosphatase, ATP diphosphohydrolase, ADP diphosphohydrolase, Adenosine 5'-diphosphatase, Adenosine 5'-triphosphatase

Many plant and animal tissues contain pyrophosphohydrolases commonly called apyrases. Apyrases catalyze the hydrolysis of a broad range of nucleoside triphosphates and diphosphates, per the following general summary reaction:^{1,2}

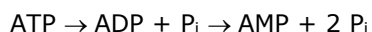

Several characteristics distinguish apyrases from other phosphohydrolases, such as:³

- High specific activity
- Broad nucleotide substrate specificity for nucleotides
- Insensitivity to specific inhibitors of P-type, F-type, and V-type ATPases

In addition, apyrases require metal cations for their activity, where the major positive effect is achieved with Ca^{+2} .

This recombinant product was cloned from *Solanum tuberosum* apyrase and expressed in *Pichia pastoris*. This isoform has similar specific activities for hydrolysis of both ATP and ADP.³ The product is lyophilized from a solution containing 50 mM potassium succinate (pH 6.5) and 100 mM NaCl.

Several publications^{4,5} and dissertations^{6,7} have cited use of product A6237 in their research protocols.

Precautions and Disclaimer

For R&D use only. Not for drug, household, or other uses. Please consult the Safety Data Sheet for information regarding hazards and safe handling practices.

Reagent

Purity: $\geq 90\%$ (SDS-PAGE)

ATPase activity: $\geq 1,000$ units/mg protein

Unit definition: One unit will liberate 1.0 μmole of inorganic phosphate from ATP or ADP per minute at pH 6.5 at 30 °C.

Preparation Instructions

This product is soluble in water, such as at 100 units/mL.⁸ It is recommended to reconstitute material in water to a concentration of 100-500 units/mL.

Storage/Stability

Store product at -20 °C. When stored at -20 °C, the enzyme retains activity for at least two years.

After reconstitution, product can be kept at 2-8 °C for up to one week. It is recommended to store the protein in working aliquots at -20 °C.⁸

References

1. Kettlun, A.M. *et al.*, *Phytochemistry*, **31(11)**, 3691-3696 (1992).
2. Komoszynski, M., and Wojtczak, A., *Biochim. Biophys. Acta*, **1310(2)**, 233-241 (1996).
3. Handa, M., and Guidotti, G., *Biochem. Biophys. Res. Commun.*, **218(3)**, 916-923 (1996).
4. Madry, C. *et al.*, *Proc. Nat. Acad. Sci. USA*, **115(7)**, E1608-E1617 (2018).
5. Zhang, X. *et al.*, *Chem. Commun. (Camb.)*, **55(98)**, 14765-14768 (2019).
6. Pezzato, Cristian, "Organization and signal regulation in complex chemical systems". Università degli Studi di Padova, Ph.D. dissertation, p. 121 (2015).
7. Aibibula, Miriayi, "Energy Metabolism of Platelets". University of Hull and University of York / Hull York Medical School, Ph.D. dissertation, p. 242 (2017).
8. Pezzato, C., and Prins, L.J., *Nat. Comm.*, **6**, 7790 (2015).

Notice

We provide information and advice to our customers on application technologies and regulatory matters to the best of our knowledge and ability, but without obligation or liability. Existing laws and regulations are to be observed in all cases by our customers. This also applies in respect to any rights of third parties. Our information and advice do not relieve our customers of their own responsibility for checking the suitability of our products for the envisaged purpose.

The information in this document is subject to change without notice and should not be construed as a commitment by the manufacturing or selling entity, or an affiliate. We assume no responsibility for any errors that may appear in this document.

Technical Assistance

Visit the tech service page at SigmaAldrich.com/techservice.

Terms and Conditions of Sale

Warranty, use restrictions, and other conditions of sale may be found at SigmaAldrich.com/terms.

Contact Information

For the location of the office nearest you, go to SigmaAldrich.com/offices.

The life science business of Merck KGaA, Darmstadt, Germany operates as MilliporeSigma in the U.S. and Canada.

MilliporeSigma, and Sigma-Aldrich are trademarks of Merck KGaA, Darmstadt, Germany or its affiliates. All other trademarks are the property of their respective owners. Detailed information on trademarks is available via publicly accessible resources.

© 2016-2022 Merck KGaA, Darmstadt, Germany and/or its affiliates. All Rights Reserved.

A6237dat Rev 10/22 DS,DT,RBG,GCY,MAM

**Millipore
Sigma**